

SWITCHING OFF X-RAY SETUP

POWER SUPPLY SETUP



STEPS:

1. Decrease the current step by step to 10 mA.
2. Decrease the voltage step by step to 10 mV.

first decrease current and then voltage .

3. Turn “OFF” the power button. (for switching off the X-ray)
4. Turn off the vacuum gauge behind the generator
5. Turn off the two vacuum pumps behind the generator

back panel



Water circulation section for cooling :

Wait for 10 minutes
for system to cool
down.

Then close the four
taps for TARGET and
TUBE only.

WARNING : the two
TMP taps need to stay
open.



Chiller :

Turn off the lowering of temperature of sample by
Press the 'OFF' switch above the chiller



TURNING OFF THE THERMOMETER :

- Turn OFF the thermometer reading the temperature of the sample.

For it turn the thermometer to PC mode and then press the OFF button.



IT'S
OVER